# Manual

## ApiPlus

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**ApiPlus is a concentrate (powder form) for the preparation of 100 ml liquid medium in which bee semen can be preserved at +17°C for 4 Wochen. ApiPlus is delivered as a powder (3.0 g) in a bottle.**

## 1. Package size

Each bottle of ApiPlus contains 3.0 g of concentrate for the preparation of 100 ml extender solution. Bi-distilled water (100 ml, Ref. 13570/1100) is required for the preparation. The water is not included in the delivery and must be ordered separately.

## 2. Application

### 2.1. Preparation of the ready to use extender

- Add 100 ml of the bidistilled water into the bottle with the powder mixture and close the bottle.
- Agitate the bottle for 5 seconds.
- Wait for 15 minutes before agitating the bottle again for another 15 seconds.
- Then wait for another 5 minutes.
- Prior to use carefully swirl the bottle once again.

### 2.2. Application for pooled ejaculates

Extend the freshly collected, pooled drone ejaculate with the extender medium depending on the semen quality and density in such a way, so that each insemination dose contains at least 4 Mio motile semen cells. We recommend a dilution ratio of 10:1 (extender to semen).

Please make absolutely sure that the semen and the extender have the same temperature.  
Avoid temperature variations during the conservation.

Slowly cool the diluted semen down to +17°C.

The diluted semen can be stored at +17°C for up to 4 weeks.

## 3. Storage, shelf life and important hints

### 3.1. Storage and shelf life

The ready made medium can be stored for 1 week at +2°C to +8°C.

For long term storage we recommend aliquot the medium and to store it at -20°C. Thus the shelf life extends to one year.

Sealed bottles can be stored dark and dry at +5°C to +10°C.

The minimum shelf life is printed on each bottle label.

### 3.2. Important hints

Keep the product out of the reach of children.

The product is not suitable for treatment or consumption by human beings or animals.

Before each use of the extender, make sure to analyse precisely, whether the extender shows the described properties, especially regarding its consistency and aspect. If the extender shows a different consistency or aspect than normal, it must not be used or only after consulting Minitüb.

### 3.3. Trouble shooting

A couple of factors are crucial for good semen conservation and insemination success, among others:

- Use of exclusively sperm friendly materials and surfaces for semen and extender.
- Avoid temperatures below +30°C for undiluted semen and temperature fluctuations.
- Avoid bacterial contamination (appropriate bee management and preparation of collection; semen collection in clean ambient, appropriate hygiene protocol in the laboratory).
- Avoid sperm toxic factors (water, UV light, dust, detergent)

## 4. Production standard and quality control

### 4.1. Raw material quality

All raw materials used in Minitube media are produced according to GMP and DIN ISO 9001:2015 norms, and are certified for meeting Ph Eur, BP or USP standards. They are tested and comply with the current international guidelines regulating the quality of pharmaceutical substances.

### 4.2. Production standard

The mixing and bottling is performed under GMP conditions in a controlled atmosphere room and documented with weighing records. Under the supervision of a team of veterinary specialists, the final product is subjected to a series of chemical and physical tests designed to ensure that Minitube media can consistently provide semen conservation with high spermatological efficacy over a wide range of climatic and process conditions.

### 4.3. Certificates

The general quality certificate and the batch certificate are available upon request.

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All rights reserved, including the translation.  
Reprint, even in extracts only with a prior written permission. No part of this manual may be reproduced or transmitted in any form or by any means electronic, mechanical, photocopying, without the prior written permission of Minitüb GmbH.  
All brand product names are trademarks or registered trademarks of the respective title holder.  
Errors and technical alterations excepted.  
Use product only in perfect condition and in compliance with the manual.  
Keep safe for future consulting (according to EN ISO 12100).  
Translated from German.

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